CAT. I

① MICROMIUM accessories

② MODEL

STERILE OVERBOOTS : WL-CBEMS-00
STERILE SLEEVE : WL-MMS-00

③ CE

Conformity to regulation EU - 2016/425

DO NOT REUSE

02 ENGLISH

03 ESPAÑOL

02 FRANÇAIS

04 DEUTSCH

ENGLISH

1. Trademark
2. Models
3. CE marking: category I PPE
4. Refer to instructions for use
5. Fabric care symbols
6. Single use

PHYSICAL PERFORMANCE OF MICROMIUM ACCESSORIES®

<i>Fabric Physical Properties</i>	<i>Test Method</i>	<i>Results</i>
Abrasion resistance	ISO 12947-2	100 cycles
Flex cracking resistance	ISO 7854 B	50 000 cycles
Trapezoidal Tear	ISO 9073-4	longitudinal direction : 44.8 N transversal direction: 15.9 N
Tensile strength	ISO 13934-1	longitudinal direction: 84 N transversal direction: 41 N
Puncture resistance	EN 863	8 N
Seam strength	ISO 13935-2	92 N

RESISTANCE TO WATER PENETRATION – INCREASING HYDROSTATIC PRESSURE

<i>Test method</i>	<i>Results</i>
ISO 811 :2018	150 cm H ₂ O

Model and area of use

White elasticated sleeve (polypropylene interior and PE exterior) with thumb loop, one size.

White overboots (polypropylene interior and PE exterior) with elasticated ankle, non-slip PVC sole, one size.

These single-use accessories are designed to provide protection against minor risks only. These accessories comply with the requirements for category I personal protective equipment as defined by Regulation (EU) 2016/425. These accessories are complementary to a category I suit. The accessories have been sterilised by gamma irradiation. They are designed in particular for work in sterile or non-sterile laboratories. The accessories are intended to protect the wearer from small splashes of cleaning materials of weak action or prolonged contact with water, and against superficial mechanical injury. They limit contamination in sensitive areas.

How to wear the accessory

How to put on the sleeve: Gently put the accessory on by inserting the arm through the folded part. Position the thumb in the elasticated loop (thumb loop).

How to put on the overboots: Gently slip the accessory on by inserting the foot through the opening, then adjust it by pulling up to the top of the calf. Tie the straps to hold and adjust the overboots.

Limit of use

The user will be the sole judge of the suitability of the protective accessories for the situation, the protection required, and the choice and combination of accessories with other equipment (gloves, respiratory protective equipment, etc.). The manufacturer cannot be held responsible for the improper or inappropriate use of the accessories. The accessories are for single use only and must be disposed after any job. Do not use if the packaging is damaged. Use the product within 3 years of the date of manufacturing (stated on the outer packaging). These accessories are for professional use only. Do not protect against biological and chemical hazards. These accessories are flammable, keep away from fire. Do not use the accessories if any defects are noticed (e.g.: holes, damaged seams, ...).

Storage and disposal

Accessories should be stored in their original packaging in a dry and away from sunlight. They must be disposed of in accordance with the regulations in force.

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MEDICOM SAS