« Work Safely
Work Together »

CAT. I

① MICROMIUM accessories

② MODEL

STERILE OVERBOOTS : WL-CBEMS-00
STERILE SLEEVE : WL-MMS-00

③ CE

Conformity to regulation EU - 2016/425

④

⑤

⑥

DO NOT REUSE

02 ENGLISH

03 ESPAÑOL

02 FRANÇAIS

04 DEUTSCH

ENGLISH

1. Brand name
2. Models
3. CE marking: category I PPE
4. Refer to instructions for use
5. Fabric care symbols
6. Single use

PHYSICAL PERFORMANCE OF MICROMIUM ACCESSORIES®

| <i>Fabric Physical Properties</i> | <i>Test Method</i> | <i>Class</i> |
|-----------------------------------|--------------------|--------------|
| Abrasion resistance | ISO 12947-2 | 2 |
| Flex cracking resistance | ISO 7854 B | 6 |
| Trapezoidal Tear MD / XD | ISO 9073-4 | 1 |
| Tensile strength MD / XD | ISO 13934-1 | 1 |
| Puncture resistance | EN 863 | 1 |
| Seam strength | ISO 13935-2 | 2 |

RESISTANCE TO WATER PENETRATION – INCREASING HYDROSTATIC PRESSURE

| <i>Test method</i> | <i>Results</i> |
|--------------------|-------------------------|
| ISO 811 :2018 | 150 cm H ₂ O |

Model and area of use

White elasticated sleeve (polypropylene interior and PE exterior) with thumb loop, one size.

White overboots (polypropylene interior and PE exterior) with elasticated ankle, non-slip PVC sole, one size.

These single-use accessories are designed to provide protection against minor risks. These accessories comply with the requirements for category I personal protective equipment as defined by Regulation (EU) 2016/425. These accessories are complementary to a category I suit. The accessories have been sterilised by gamma irradiation. They are designed in particular for work in sterile or non-sterile laboratories. They are designed to protect the wearer from small splashes, and to limit contamination in sensitive areas.

How to wear the accessory

How to put on the sleeve: Gently put the accessory on by inserting the arm through the folded part. Position the thumb in the elasticated loop (thumb loop).

How to put on the overboots: Gently slip the accessory on by inserting the foot through the opening, then adjust it by pulling up to the top of the calf. Tie the straps to hold and adjust the overboots.

Limit of use

The user will be the sole judge of the suitability of the protective accessories for the situation, the protection required, and the choice and combination of accessories with other equipment (gloves, respiratory protective equipment, etc.). The manufacturer cannot be held responsible for the improper or inappropriate use of the accessories. The accessories are for single use only and should be disposed of after use. Do not use if the packaging is damaged. Use the product within 3 years of the date of manufacture (stated on the outer packaging). These accessories are for professional use only. Do not protect against biological and chemical hazards.

Storage and disposal

Accessories should be stored in their original packaging in a dry and dark place. They must be disposed of in accordance with the regulations in force. Restrictions on their disposal depend solely on the pollutants that may have come into contact with the garment during use.

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