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## ENGLISH

1. Brand name
2. Model identification
3. CE marking : Category III PPE
4. European Protection Standards
5. Size and body measurements
6. Refer to instructions for use
7. Fabric care symbols
8. Single use

## PHYSICAL PERFORMANCE OF WEEPRO LABO®

Fabric Physical Properties	Test Method	Class
Resistance to ignition	EN 13274-4	PASS
Abrasion resistance	ISO 12947-2	3
Flex cracking resistance	ISO 7854 B	6
Impact resistance	ISO 9073-4	1
Tensile strength MD7/XD	ISO 13934-1	1
Truncture resistance	EN 863	1
Seam strength	ISO 13935-2	3

## RESISTANCE TO CHEMICAL LIQUID EN ISO 6530

Fabric Chemical Properties	Penetration	Repellency
Sulphuric Acid 30 %	CLASS 3	CLASS 3
Sodium Hydroxide 10 %	CLASS 3	CLASS 3
Butan-1-ol	CLASS 3	CLASS 3
o-xylene	CLASS 3	CLASS 2

## WHOLE SUIT TEST PERFORMANCE

Dry particles Type 5 Method as defined by EN ISO 13982-1:2004 + A1:2010 | PASS  
Light splashes Type 6 Method as defined by EN 13034 : 2005 +A1: 2009 | PASS

## Model and typical area of use

White coveralls with elasticated hood, wrists, ankles and back, adhesive flap over the zipper. Coveralls with integrated boots. Coveralls are designed to protect against chemical risks (splashes – light sprays – airborne solid particulates).

## How to wear protective clothing

Remove the coverall from its packaging, open the center zipper and put it on the garment, taking care not to damage it. Close the zipper completely and cover it with the adhesive flap, without creasing. If there is a risk of airborne solid particles, it is advisable to tape additionally protective gloves with ends of the sleeves.

## Limitations of use

Do not use the coverall if any defects are noticed (e.g.: holes, damaged seams and fastenings, ...). The coverall meets the requirements:  $L_{min} 62/90 \leq 30\%$  -  $L_{s, 8/10} \leq 15\%$  % Type 6 chemical protective suits have been tested to the whole suit test. Wear for long periods of time can cause heat stress. Heat stress and discomfort can be reduced or eliminated by using appropriate undergarments or suitable ventilation equipment. Coveralls are for single use only and must be disposed after each use. The user is the sole judge of the adequacy between his risk analysis and the type of protection required, including the additional equipment necessary for a complete protection of the body. If necessary, use additional PPE with the same level of protection (such as gloves, breathing apparatus, etc.) in order to provide full body protection. WeeSafe® cannot accept responsibility for any improper use or disposal of garments produced by them. These garments are flammable, keep away from fire. Use the product within 3 years of the date of manufacturing.

## Storage and disposal

The coveralls must be stored in their original packaging in a dark, dry place. They must be disposed of in compliance with current regulations.

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